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liquitec stands for stringent standards of quality, reliability, engineering and process design

#### About liquitec AG

Safety assured by more than 25 years of experience and know-how:

- Aseptic design is our standard
- Qualifiable for use in your cGMP plants
- Exceptional cleanability reduces CIP costs
- Hermetically sealed for absolute sterility
- Scalability enables down and up scaling

Typical applications for our products include:

#### Bio-Technology

- Bioreactors for bacteria and cell cultures
- Upstream and downstream process vessels
- Vaccine production

#### Pharma-Technology

- active pharmaceutical ingredient
- Infusion and injection solutions

#### Food-Technology

- Fruit juice production
- Breweries
- Dietary supplements
- Sterile vessels

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Achieve maximum flexibility and economy with our modular system bio-m<sup>®</sup>, bio-plus<sup>®</sup>, bio-t<sup>®</sup> & bio-pure<sup>®</sup>

Our modular system bio-m<sup>®</sup> enables our customers to achieve maximum flexibility and cost-effectiveness in their process designs, plant construction and plant maintenance concepts. The latest technological advances and uncompromising sterile designs are key differentiators of all our bio-m<sup>®</sup> products.

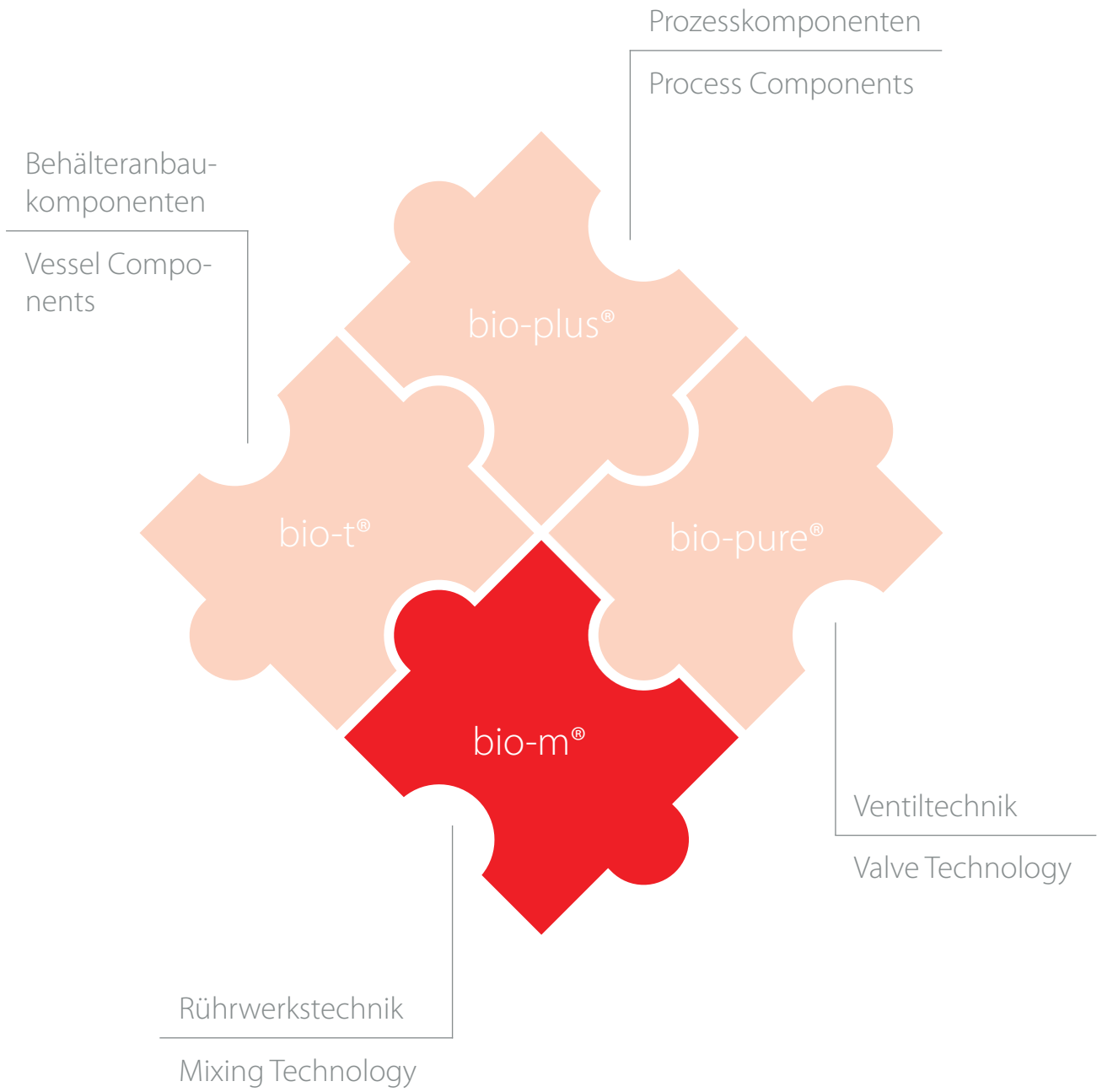
bio-plus<sup>®</sup> is the name for our program of process components like pipe unions, strainers, sight glasses or non-return valves.

Our bio-t<sup>®</sup> component system offers a complete, high-quality product range of vessel components.

bio-pure<sup>®</sup> sterile valve technology combines the advantages of the classic diaphragm valve with numerous, partly patented, innovative product characteristics. bio-pure<sup>®</sup> is characterized by strict sterile design, compact construction, flexibility in application and shortest maintenance times!

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
Sterile technology within a System

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
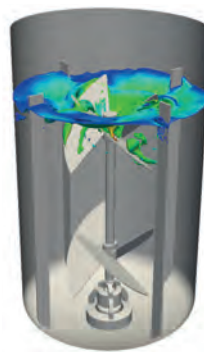
## Fundamentals of Mixing Technology

Mixing is one of the most important basic operations of applied processing technology. In nearly all processes in the chemical, pharmaceutical, biological and food industries, as well as in raw material and water preparation, mixing is used to ensure homogenization or heat and mass transfer processes in liquid systems.

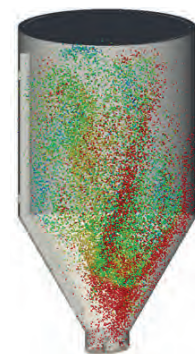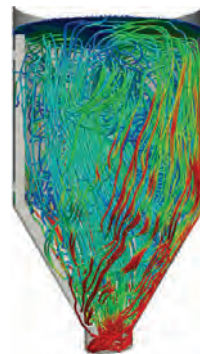The job of mixing is to create a three-dimensional flow pattern in the container with high pulse current densities, as well as achieving the largest possible boundary surface in multiphase systems. Optimum mixer design requires the latest scientific methods (e.g. CFD simulations) as well as practical experience.

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## Mixing Tasks

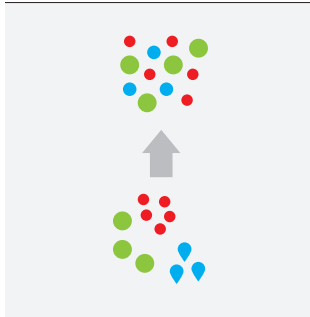
liquitec designs machines to suit the required process functions. The typical functions listed below can be performed either separately or in combination.

- Homogenizing
- Suspension
- Heat transfer
- Gas flushing
- Dispersing

With such a large number of critical mixing functions to perform, it is important to seek competent advice on optimal and safe process control. Our mixer designs are matched to your application based on your process description and specifications.


On request, we can also analyse your existing mixing processes and provide suggestions regarding any potential for optimization that may exist.

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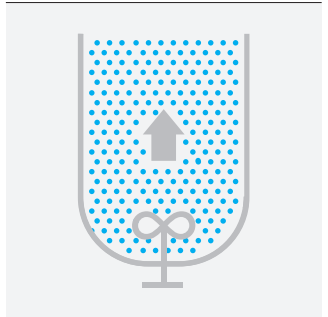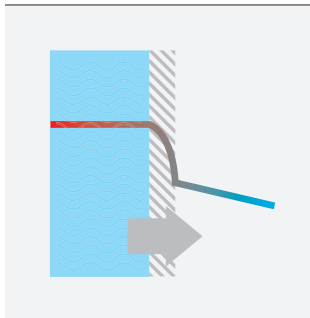
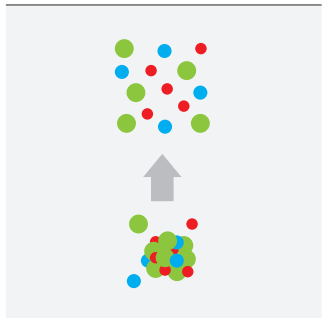
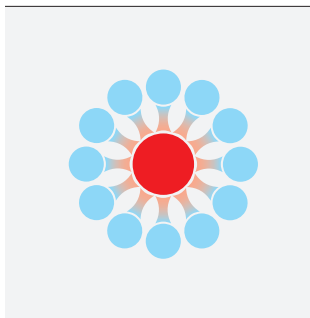
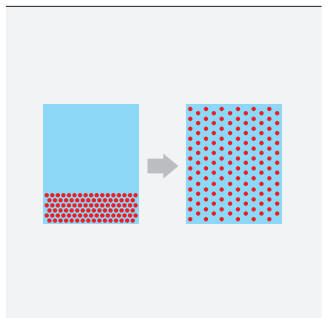
## Directives and Certificates

### liquitec Quality

The continuous development and improvement of our processes and steady investment in new production technologies help to guarantee the high quality for all liquitec products. All production steps required in the manufacturing of liquitec products, are completely documented and can be traced at any time. Our process oriented internal quality management helps to provide our customers with the best products and services.

Regarding quality, liquitec also understands the knowledge and application of all relevant guidelines. These include for example ATEX, EC Machinery Directive, ASME BPE or proof of cleaning ability according to the EHEDG procedure. We only use materials with corresponding material certificates, be it according to DIN / EN 10204 or elastomers with FDA / USP Class VI / ADI-free certificates. This ensures that our customers receive high-quality and qualified agitators for their applications.

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Made of 1.4435 / AISI 316L stainless steel, Hastelloy or other specialized alloys

#### **Materials**

All liquitec product-contact parts are made of 1.4435 / AISI 316L stainless steel, usually to directive BN2, which can be certified and traced back to its base at any time. Thus, the controlled ferrite content of the materials is guaranteed at all times.

All our product-contact surfaces comply with normal industry standards and are manufactured to  $Ra < 0.6$  microns, which can be decreased to  $Ra < 0.2$  microns as an option.

liquitec products are treated in our own in-house electro-polishing facility and are packaged and delivered to our customers with chemically pure surfaces.

Alternative material options such as Hastelloy or other specialized alloys are available on request for all liquitec products.

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FDA  
USP Class VI (121°C)  
REACH / RoHS  
Sanitary 3A  
ADI-Free

**Certificates for  
elastomer sealings**

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**Business Areas**

Pharma  
Chemistry  
Biotech  
Food  
Machine and plant construction  
Special machines

Grosse Spaltmasse  
Large clearance

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### Certification of components

Material certification acc. EN 10204  
Ferrite content  
Surface quality  
Electro-polishing certification  
Endoscopy  
Weld certification  
Dye penetration / X-ray examination  
Pressure test protocol

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Innovative use of construction materials

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### Advantages

Improved energy efficiency  
Reduced cleaning costs  
Less cleaning time  
Low maintenance costs

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## Safe, contamination-free processing

Our magnetic mixers represent the culmination of 25 years of experience and continual development. Accordingly, they are the current state of the art. Their advantages for your sterile process are:

### Advantages for your aseptic processing

- Open, easily-cleaned design
- Large gap between the magnetic housing and magnetic impeller for better cleaning and greater process safety
- Advanced ceramic slide bearings bushings for high reliability and long service life
- Dissimilar materials reduce friction
- Inner slide bearing made of impact-resistant zirconium oxide
- External slide bearing made of silicon carbide: can be replaced on site if necessary
- Standard material 1.4435/AISI 316L, surface finish  $Ra \leq 0.6 \mu m$ , electrochemically polished. Other materials, higher-grade surface finishes or defined ferrite contents upon request
- Speed sensor, an optional feature, for non-contact speed monitoring of the magnetic rotor
- Mixer blade shapes determined by comparative computational fluid dynamics (CFD) tests: short mixing times, low residual miscible volume, high efficiency with optimal ease of cleaning

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Bottom Magnetic Agitator with low shear force impact

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### Standard Sizes with bio-m® standard impeller

The bio-m® standard agitator is characterized by an extremely wide range of applications. It can be used for various stirring tasks in containers from 2 to 80'000 liters.

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### Standard Sizes with bio-m® high-efficiency impeller

The high-efficiency agitator element is characterized by an extremely low shearing force impact on the product. Its versatility suits it to various process applications!

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Bottom Magnetic Agitator with medium shear force impact

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### Standard Sizes with bio-m® Rotek impeller

The Rotek impeller is a perfect option for many upstream processes. The agitator generates medium shear forces and generates a high energy input. It is ideal for dissolving difficultly soluble powders in liquids.

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Bottom Magnetic Agitator with high shear force impact

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### Standard sizes with bio-m® rotor-stator impeller

The bio-m® rotor-stator system provides high shear forces with optimal product solvability. Magnetically driven, the system always guarantees absolute sterility. It is ideal for dissolving difficultly soluble powders in liquids.

Typ BMA Type BMA	Rührer Ø [mm] Agitator Ø [mm]	P [kW] P [kW]
30 RS	52	0.18
100 RS	85	0.37
550 RS	110	1.50
2500 RS	165	5.50

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BMAF-Bottom Magnetic Agitator for installation in Fermenters

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Bottom-mounted magnetic mixers used in bioreactors allow long, contamination-free process cycles, especially when it comes to cell culture fermentation.

Given the stable design of the slide bearing, our BMAF mixers allow long, free-flying shafts, and are presently used in bioreactors of up to 15,000 L process volume. Impellers are chosen to suit the process and customer specifications. Common designs include segmented, pitched blade and Rushton impellers.

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## bio-m® Top-mounted Magnetic Agitators

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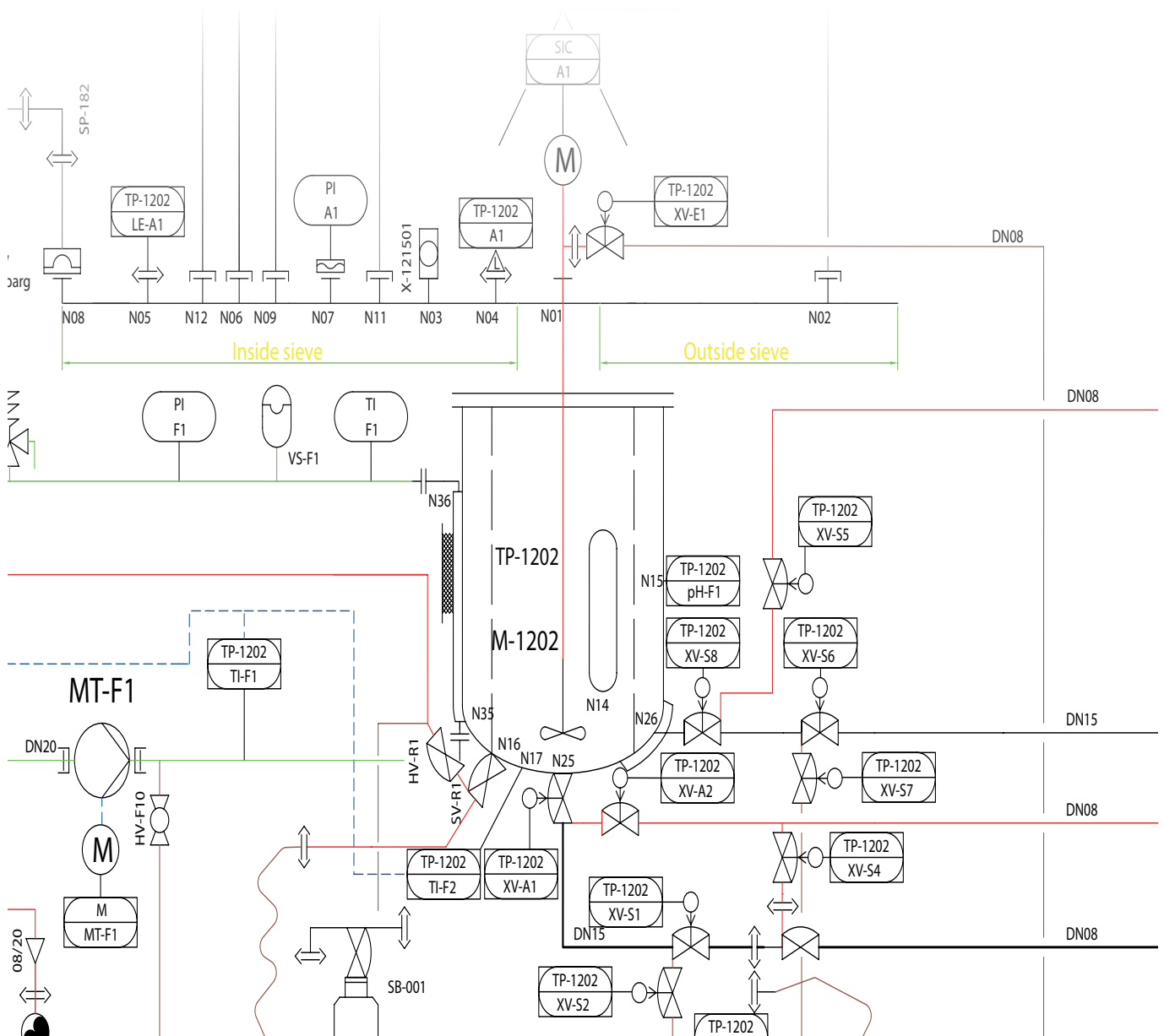
### Hermetically-sealed design assures aseptic processing

The two TMAi and TMAe model series are liquetec's solution for applications that require a magnetic mixer mounted on the vessel lid.

#### Advantages for aseptic processing


- Magnetic coupling eliminates mechanical seal
- No complicated supply systems requiring qualification, low operating costs
- Can be mounted onto an existing vessel flange if, e.g., an existing mechanical seal mixer is to be replaced
- All impeller types possible

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bio-m® TMAi Top-mounted Magnetic  
Agitator with internal Bearing


Typ TMAi Type TMAi	Drehmoment [Ncm] Torque [Ncm]
30	30
100	100
300	300
550	550

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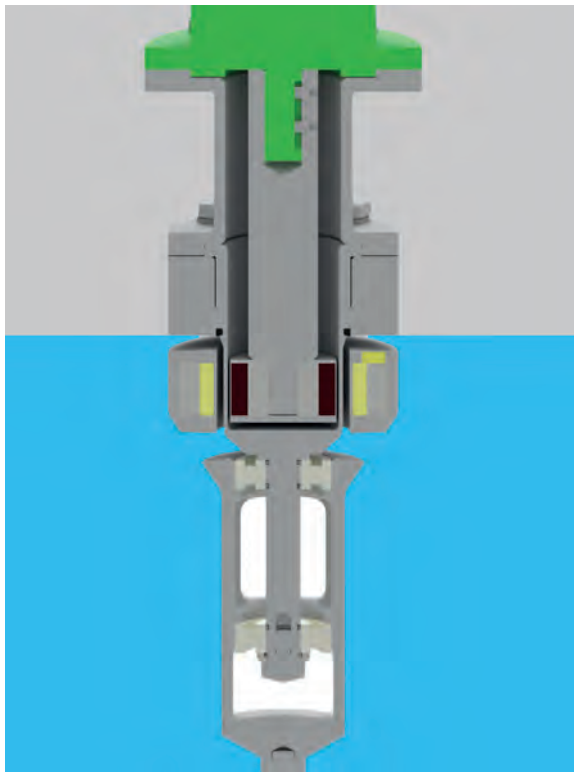bio-m® TMAe Top-mounted Magnetic Agitator  
with external Bearing

Typ TMAe Type TMAe	Drehmoment [Ncm] Torque [Ncm]
300	300
550	550
850	850
1800	1800
2500	2500

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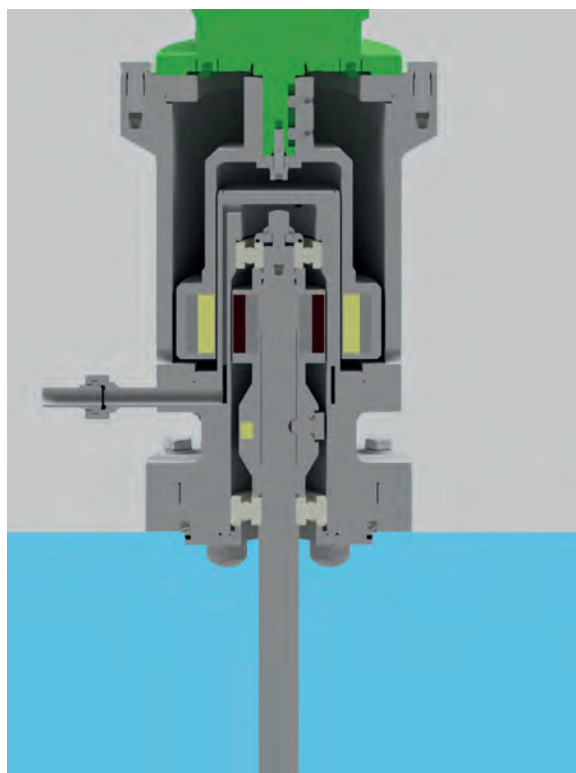### bio-m® BSA Bottom-mounted Agitator with Mechanical Seal

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These mixers with standard shaft diameters from 40 mm to 100 mm are mounted from below in the base of the vessel.

#### Vessel Connection

Exclusively dead-leg-free insertion flange, fully drainable. Drive / bearing: appropriately-dimensioned geared motors with hollow shafts and bearings integrated into the mechanical seal ensure smooth and quiet running of the mixer shaft.

Optional lowering device: allows installation of the mechanical seal while the drive is still mounted. Especially in bottom-mounted mixers, a lowering system can be a great help when installing or removing the mechanical seal.

#### Application

BSA type mixers are often used in bioreactors. The design of the impellers and speeds are defined specifically for the respective process (e.g. cell cultures or bacterial fermentation).

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bio-m® TSA Top-mounted Agitator with Mechanical Seal

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These mixers with standard shaft diameters from 25 mm to 100 mm are top-mounted on the lid of the vessel.

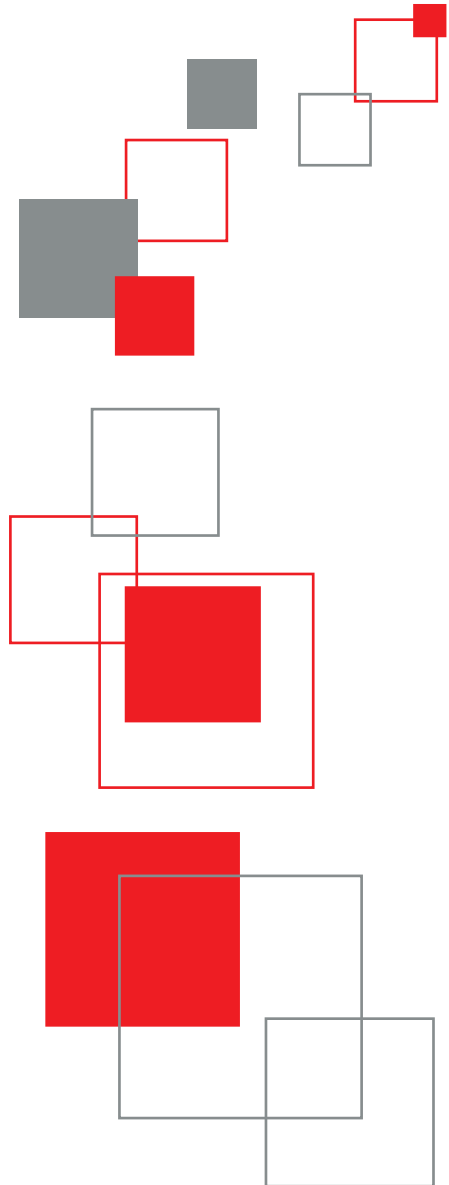
### **Container Connection**

Typically made with a dead-leg-free insertion flange. Mounting flanges according to DIN 28141 are a possible alternative.

### **Drive / Bearing**

Appropriately-dimensioned geared motors with hollow shafts and bearings integrated into the mechanical seal ensure smooth and quiet running of the mixer shaft.

 liquitec  

liquitec AG  
Industrie Neuhof 54  
CH-3422 Kirchberg  
T +41 55 450 83 00  
F +41 55 450 83 01  
[info@liquitec.ch](mailto:info@liquitec.ch)  
[www.liquitec.ch](http://www.liquitec.ch)